Digital Lock-In Amplifiers

SR850 — DSP lock-in amplifier with graphical display



• 1 mHz to 102.4 kHz frequency range

- >100 dB dynamic reserve
- 0.001 degree phase resolution
- Time constants from 10 µs to 30 ks (up to 24 dB/oct rolloff)
- Auto-gain, -phase, -reserve and -offset
- Data logging (up to 65k samples)
- Smoothing, curve fitting & statistics
- GPIB, RS-232 and 3.5" disk drive

SR850 DSP Lock-In Amplifier

The SR850 is a digital lock-in amplifier based on an innovative DSP (Digital Signal Processing) architecture. The SR850 boasts a number of significant performance advantages over traditional lock-in amplifiers—higher dynamic reserve, lower drift, lower distortion, and dramatically higher phase resolution. In addition, the CRT display and 65,536 point memory make it possible to display and process data in a variety of formats unavailable with conventional lock-ins.

Digital Precision

At the input of the SR850 is a precision 18-bit A/D converter which digitizes the input signal at 256 kHz. The A/D converter, together with a high-speed DSP chip, replace the analog demodulator (mixer), low-pass filters and DC amplifiers found in conventional lock-ins. Instead of using analog components, the SR850 is implemented by a series of precise mathematical calculations which eliminate the drift, offset, non-linearity and aging inherent in analog components. The same DSP chip digitally synthesizes the reference oscillator, providing a source with less than -80 dBc distortion, 100 mHz frequency resolution and 2 mV of amplitude resolution.


Digital Flexibility

The SR850 has a 7" CRT display which supports a large selection of display options. Data can be viewed numerically or graphically in bar graph, polar plot, and strip chart formats. With 65,536 points of memory and data acquisition rates up to

www.gmp.ch

GMP SAMain office: Avenue des Baumettes 17CH -1020 RenensTél. 021 633 21 21Fax. 021 633 21 29info@gmp.chGMP SABüro Zürich: Dübendorfstrasse 11aCH-8117 FällandenTel. 044 825 34 00Fax. 044 825 34 01info@gmp.ch

512 Hz, you are able to see exactly how your data changes in time—not just what the current output value is. After data has been acquired, the SR850 offers a variety of data reduction options, such as Savitsky-Golay smoothing, curve-fitting, and statistical analysis. A built-in 3.5" disk drive, along with standard RS-232 and GPIB interfaces, makes it easy to transfer data to your computer.

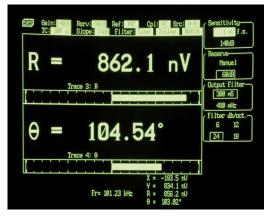

Large numeric readout with bar graph

Input Channel

The SR850 has a differential input with 6 nV/ $\sqrt{\text{Hz}}$ input noise. The input impedance is 10 M Ω , and minimum full-scale input voltage sensitivity is 2 nV. The input can also be configured for current measurements with selectable current gains of 10⁶ and 10⁸ V/A. A line filter (50 Hz or 60 Hz) and a 2× line filter (100 Hz or 120 Hz) are provided to eliminate line related interference. However, unlike conventional lock-in amplifiers, no tracking band-pass filter is needed at the input of the SR850. This filter is used by conventional lock-ins to increase dynamic reserve. Unfortunately, band-pass filters also introduce noise, amplitude and phase error, and drift. The DSP based design of the SR850 has such inherently large dynamic reserve that no tracking band-pass filter is needed.

Reference Channel

The reference source for the SR850 can be an externally applied sine wave or square wave, or its own digitally synthesized reference source. Because the internal reference source is synthesized from the same digital signal that is used to multiply the input, there is virtually no reference phase noise when using the internal reference. The internal reference can operate at a fixed frequency or can be swept linearly or logarithmically over the entire operating range of 1 mHz to 102.4 kHz. Harmonic detection can be performed at any integer harmonic of the reference frequency—not just the first few harmonics.

The DSP approach also offers considerable advantages when working with an external reference. The time to acquire an external reference is only 2 cycles + 5 ms (or 40 ms, whichever is greater)—about ten times faster than conventional lock-ins.

Because the SR850 uses a digital phase-shifting technique rather than analog phase-shifters, the reference phase can be adjusted with one millidegree resolution. In addition, the X and Y outputs are orthogonal to within one millidegree.

Outputs and Time Constants

The output time constants on the SR850 are implemented digitally. Low-pass-filter rolloffs of 6, 12, 18 and 24 dB/octave are available, with time constants ranging from 10 μ s to 30 ks. Below 200 Hz, the SR850 can perform synchronous filtering. Synchronous filters notch out multiples of the reference frequency—an especially useful feature at low frequencies where the proximity of the 2f component would otherwise require a long time constant for effective filtering. The SR850 makes working at low frequencies a far less time consuming task.


High Dynamic Reserve

The dynamic reserve of a lock-in amplifier at a given fullscale input voltage is the ratio (in dB) of the largest interfering signal to the full-scale input voltage. The largest interfering signal is defined as the amplitude of the largest signal at any frequency that can be applied to the input before the lock-in cannot measure a signal with its specified accuracy.

The SR850 has the highest dynamic reserve (>100 dB) of any lock-in available. In conventional lock-in amplifiers, dynamic reserve is increased at the expense of stability. Because of the digital nature of the filtering and gain process in the SR850, the ultra-high dynamic reserve is obtained without any sacrifice in stability or accuracy. In addition, the SR850's high dynamic reserve is obtained without the use of analog band-pass filters, eliminating the noise and error that such filters introduce.

Traces and Displays

Data acquired by the SR850 is stored in up to four userdefined traces. Each trace can be configured as $(A \times B)/C$, where A, B and C are selected from X, Y, R, θ , noise, frequency or any of the four rear-panel auxiliary inputs. Common

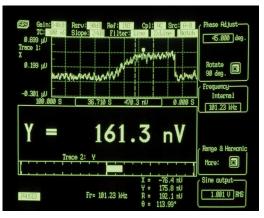

Graphical, numerical, and bar-graph display

www.gmp.ch

GMP SAMain office: Avenue des Baumettes 17GMP SABüro Zürich: Dübendorfstrasse 11a

CH -1020 Renens CH-8117 Fällanden

Tel. 044 825 34 00

Tél. 021 633 21 21

Fax. 021 633 21 29 info@gmp.ch Fax. 044 825 34 01 info@gmp.ch operations, such as ratioing, can be performed in real time by defining an appropriate trace. Trace values can be displayed as a bar graph with an associated large numerical display, or as a strip chart showing the trace values as a function of time. Additionally, you can display polar plots showing the phasor



Polar plot display

formed by the in-phase and quadrature components of the signal. All displays can be easily scaled from the front panel or over the computer interfaces, and an auto-scale feature is available to quickly optimize the display. The screen can be configured as a single large display, or as two horizontally-split displays.

Convenient Auto Measurements

Common measurement parameters are available as single-key "auto" functions. The gain, phase, dynamic reserve and display scaling can all be set with a single key press. For many measurements, the instrument can be completely configured simply by using the auto functions.

Auxiliary A/Ds and D/As

Four rear-panel A/D inputs allow you to measure external signals with millivolt resolution. The measured values can be incorporated into one of the SR850's trace definitions, displayed on the front panel, or read via either computer interface. Four D/A outputs can provide either fixed output voltages or a voltage level which scans synchronously with the SR850's frequency scans. Both the A/D inputs and the D/A outputs have a ± 10 V range.

Analysis Features

The SR850's performance doesn't stop once data has been acquired—a full set of data processing features is also included. Multiple-range Savitsky-Golay smoothing can be applied to any of the trace arrays, and statistical information (mean, variance, sum) can be calculated for a selected trace region. A curve fitting routine calculates best fits to lines, exponential curves, and Gaussian curves for any portion of your data. And a trace "calculator" lets you perform a variety of simple arithmetic and trigonometric operations on trace data.

Interfaces and Hardcopies

The SR850 comes standard with RS-232 and GPIB interfaces. All instrument functions can be queried and controlled via the interfaces. For convenient debugging, characters received and sent via the interfaces can be viewed on the front panel. Several hardcopy options are available on the SR850. Screens can be dumped to a dot-matrix or LaserJet compatible printer through the standard Centronics printer interface. Displays can also be plotted on any HP-GL compatible plotter via GPIB or RS-232.

Ordering	Information
SR850	DSP dual phase lock-in
	amplifier (w/ rack mount)
O850H	Carrying handle kit
SR550	Voltage preamplifier
	$(100 \text{ M}\Omega, 3.6 \text{ nV}/\sqrt{\text{Hz}})$
SR552	Voltage preamplifier
	$(100 \text{ k}\Omega, 1.4 \text{ nV}/\sqrt{\text{Hz}})$
SR554	Transformer preamplifier
	(0.091 nV/\Hz)
SR540	Optical chopper



www.gmp.ch

GMP SAMain office: Avenue des Baumettes 17GMP SABüro Zürich: Dübendorfstrasse 11a

 CH -1020 Renens
 Tél. 021 633 21 21
 Fax. 021 633 21 29
 info@gmp.ch

 CH-8117 Fällanden
 Tel. 044 825 34 00
 Fax. 044 825 34 01
 info@gmp.ch

SR850 Specifications

Signal Channel

Voltage inputs Sensitivity Current input Input impedance Voltage input Current input Gain accuracy Noise (typ.)

Line filters CMRR

Dynamic reserve

Reference Channel

Frequency range Reference input Input impedance Phase resolution Absolute phase error Relative phase error Orthogonality Phase noise Int. reference Ext. reference Phase drift

Harmonic detection Acquisition time

Demodulator

Stability Digital outputs Analog outputs Harmonic rejection Offset/Expand Time constants

Internal Oscillator

Range Accuracy Resolution

Distortion

Amplitude

Output impedance Amplitude accuracy Amplitude stability Outputs

Sweeps

Single-ended or differential 2 nV to 1 V 10^6 or 10^8 V/A $10 \text{ M}\Omega + 25 \text{ pF}$, AC or DC coupled 1 k Ω to virtual ground

±1 % (±0.2 % typ.) $6 \text{ nV}/\sqrt{\text{Hz}}$ at 1 kHz $0.13 \text{ pA}/\sqrt{\text{Hz}}$ at 1 kHz (10⁶ V/A) $0.013 \text{ pA}/\sqrt{\text{Hz}}$ at 100 Hz (10⁸ V/A) 50/60 Hz and 100/120 Hz (Q=5) 100 dB at 10 kHz, decreasing by 6 dB/oct above 10 kHz >100 dB (without prefilters)

0.001 Hz to 102.4 kHz TTL or sine (400 mVpp min.) 1 MΩ, 25 pF 0.001° <1° <0.001° $90^{\circ} \pm 0.001^{\circ}$ <0.0001° rms at 1 kHz 0.005° rms at 1 kHz, 100 ms, 12 dB/oct

<0.01°/°C below 10 kHz, <0.1°/°C, 10 kHz to 100 kHz 2F, 3F, ... nF to 102.4 kHz (2 cycles + 5 ms) or 40 ms,whichever is greater

no drift <5 ppm/°C for all dynamic reserves –90 dB ± 100 % offset, expand up to 256× 10 µs to 30 ks (6, 12, 18, 24 dB/oct rolloff) Synch. filtering available below 200 Hz.

1 mHz to 102.4 kHz $25 \text{ ppm} + 30 \mu\text{Hz}$ 0.01 % or 0.1 mHz (whichever is greater) -80 dBc (f < 10 kHz)-70 dBc (f > 10 kHz) at 1 Vrms 0.004 to 5 Vrms into 10 k Ω (2 mV resolution) 50 Ω 1 % 50 ppm/°C Sine and TTL (both can be phaselocked to an external reference) Linear and log

Inputs and Outputs

inputs and outputs	
Interfaces	IEEE-488.2, RS-232 and Centronics interfaces standard. All instrument functions can be controlled and read though the interfaces
X, Y outputs CH1 output CH2 output Aux. A/D inputs Aux. D/A outputs	though the interfaces. ± 10 V, updated at 256 ksamples/s ± 10 V output of X, R or Trace 1 to 4 ± 10 V output of Y, θ or Trace 1 to 4 4 BNC inputs, 1 mV res., ± 10 V 4 BNC outputs, 1 mV resolution, ± 10 V (fixed or swept amplitude)
Sine out TTL out Trigger In	Internal oscillator analog output Internal oscillator TTL output TTL signal starts internal oscillator sweep or triggers instrument
Remote pre-amp	data taking (rates to 512 Hz). Provides power to the optional SR550, SR552 and SR554 preamps
Displays	
Screen format Displayed quantities	Single or dual display Each display shows one trace. Traces are defined as $A \times B/C$ or $A \times B/C^2$ where A, B, C are selected from X, Y, R, θ , X-noise, Y-noise, R-noise,
Display types	Aux 1 to 4 or frequency. Large numeric readout, bar graph, polar plot and strip chart
Data buffer	64k data points. Buffer is configured as a single trace with 64k points, two traces with 32k points each, or four traces with 16k points each.
Sample rate	0.0625 to 512 Hz, external to 512 Hz
Analysis Functions	
Smoothing Curve fitting Calculator	5, 9, 17, 21, 25 pt. (Savitsky-Golay) Linear, exponential or Gaussian Arithmetic, trigonometric and logarithmic calculations
Statistics	Mean and standard deviation
General	
Hardcopy	Screen dumps to dot-matrix or LaserJet printers. Plots to HP-GL plotters (RS-232 or GPIB).
Disk drive	3.5" MS-DOS compatible format, 1.44 Mbyte. Storage of data and instrument setups (binary or ASCII). Screens can be saved as PCX files.
Power	60 W, 100/120/220/240 VAC, 50/60 Hz
Dimensions	17" × 6.25" × 19.5" (WHD)
Weight	40 lbs.
Warranty	One year parts and labor on defects in materials and workmanship

www.gmp.ch

GMP SA Main office: Avenue des Baumettes 17 GMP SA Büro Zürich: Dübendorfstrasse 11a

CH -1020 Renens CH-8117 Fällanden

Tel. 044 825 34 00

Tél. 021 633 21 21 Fax. 021 633 21 29 Fax. 044 825 34 01

info@gmp.ch info@gmp.ch

SR850 DSP Lock-In Amplifier

About DSP Lock-In Amplifiers

DSP lock-in amplifiers differ considerably from their analog counterparts as a glance at the block diagram below will quickly reveal. Although the front-end of both types of lock-ins contain a low-noise AC amplifier and line filters, the similarity ends there. In the SR850, the signal is filtered with a 108 kHz, 9th order, elliptical anti-aliasing filter. This filtering is crucial to ensure that the signal can be digitized by the 256 kHz, 18-bit A/D converter with no aliasing.


The A/D converter passes the digitized signal to the DSP chip. The DSP chip synthesizes a 24-bit digital reference sine wave at the reference frequency. In the internal reference case, the reference is synthesized from a high-accuracy crystal oscillator. In the external reference case, a phase-locked loop locked to the external reference serves as a source for the reference signal. The reference is multiplied by the signal in the DSP chip, which is capable of performing 16 million 24-bit \times 24-bit multiplies and additions each second. After multiplication, up to four stages of digital low-pass filtering are applied to generate time constants from 10 µs to 30 ks, with filter rolloffs of 6, 12, 18, and 24 dB/octave.



The resulting X and Y (in-phase and quadrature) signals are used to digitally calculate the values of R and θ . The results are sent to the main system processor for display on the CRT, and passed through an 18-bit digital-to-analog converter to generate the front-panel outputs.

The same digital reference that was used to multiply the signal is converted by another 18-bit D/A converter, and is used as the SR850's internal oscillator. Thus, the internal oscillator output is actually the same signal as the reference, and there is virtually no reference phase noise when using it. Phase shifting the reference is also simple in this model, since only digital calculations are involved.

It is this digital architecture which makes possible the performance advantages found in the SR850. Comparing the diagram below with a block diagram of a conventional lock-in shows that many of the most troublesome components—noisy input prefilters, nonlinear demodulators, inaccurate analog filters, and drift-prone, high-gain DC amplifiers—have all been removed, along with their performance penalties. The resulting instrument comes closest to implementing the theoretical model of lock-in amplification.

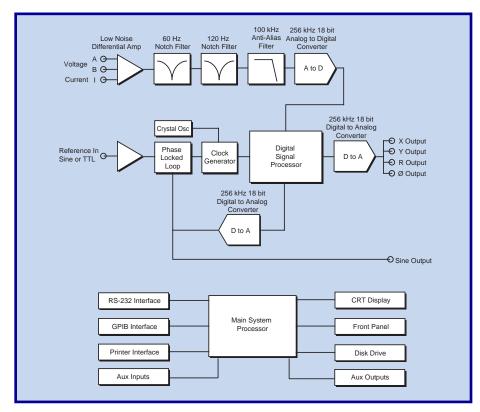

DSP Lock-In Amplifier Block Diagram

www.gmp.ch

GMP SA Main office: Avenue des Baumettes 17GMP SA Büro Zürich: Dübendorfstrasse 11a

CH -1020 RenensTél. 021 633 21 21CH-8117 FällandenTel. 044 825 34 00

 Fax. 021 633 21 29
 info@gmp.ch

 Fax. 044 825 34 01
 info@gmp.ch